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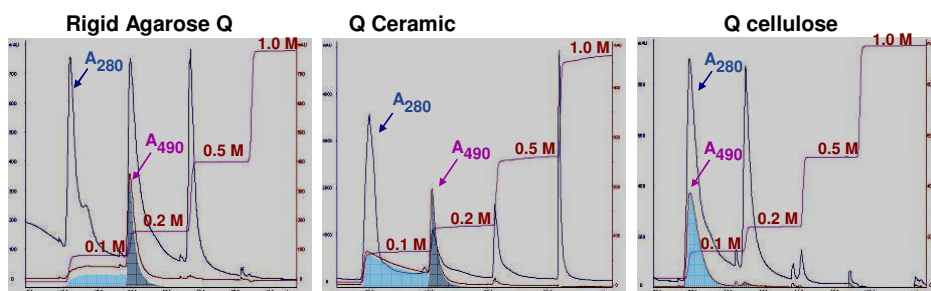
Purification of rGFP from E.coli lysate on anion exchangers

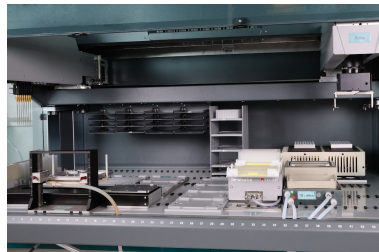
50 mM Tris-HCl pH 8.5; step elution at 0.1, 0.2, 0.5 and 1M NaCl; 0.5 cm I.D. x 5 cm at 300 cm/h (1 min RT)

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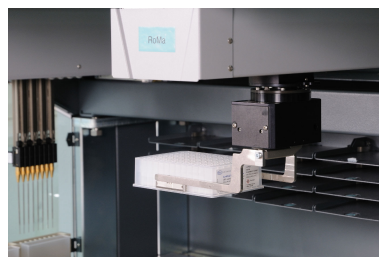
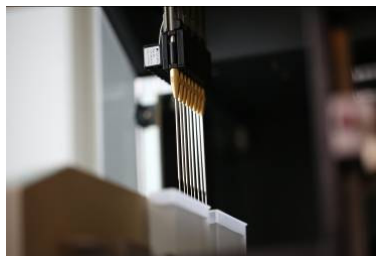
PALL Life Sciences

2013: Pall ScreenExpertSM Services Laboratory Cergy (Paris), France

7

PALL Life Sciences

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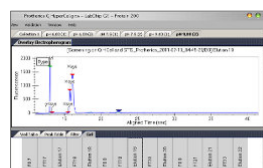
SELDI-Mass Spectrometry

Automated microfluidic
electrophoresis (LabChip®)

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

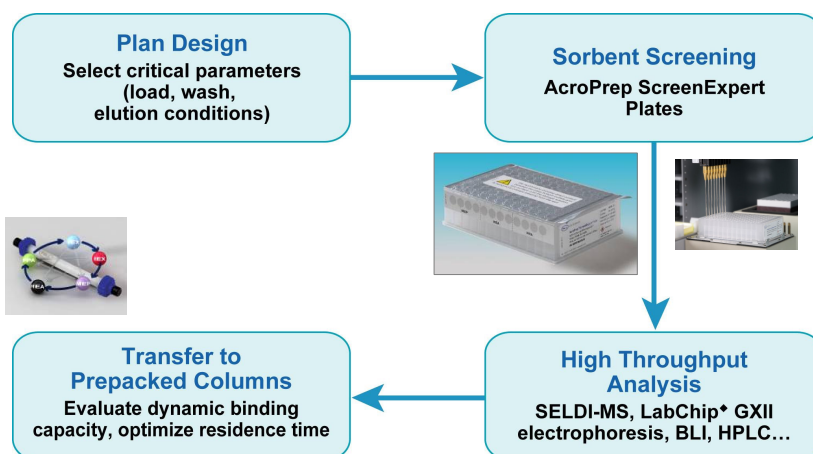
11

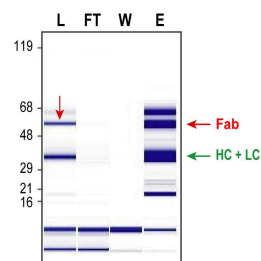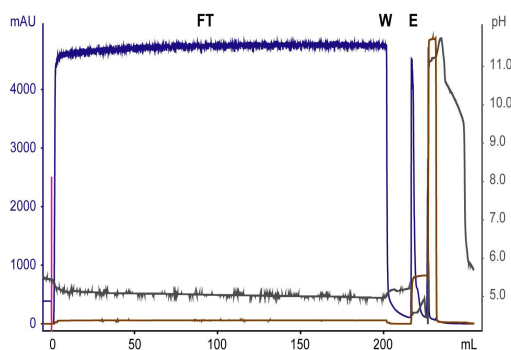Capture step by Cation exchange: CM Ceramic HyperD® 1 mL PRC Column

Capture step performance

- Fab recovery = 96%
- Concentration = 20 times
- Purity = 40%

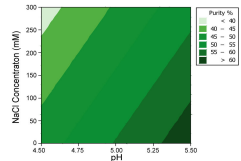
Equilibration: 50 mM Na acetate pH 5.0
Loading: 200 mL feedstock 1/2 at pH 5.0
Washing: 50 mM Na acetate pH 5.0
Elution: 50 mM Na acetate pH 5.0 + 0.3M NaCl
Flow rate: 150 cm/h – Residence Time = 3 min

12



Intermediate step: Screening on AcroPrep 96 plates (Mixed Mode on MEP HyperCel)

Fab purity

Fab recovery

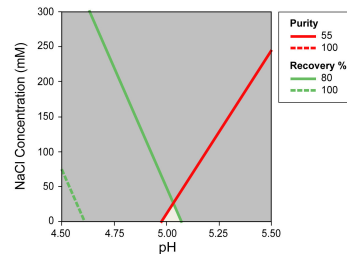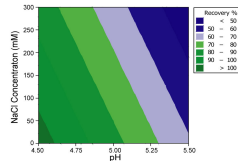
Screening of 3 pH (pH 4.5, 5.0, 5.5) and
3 NaCl conditions (0 mM, 150 mM, 300 mM)



Optimal conditions on MEP HyperCel (FT mode):
Binding at pH 5.0 to 5.2, without addition of NaCl

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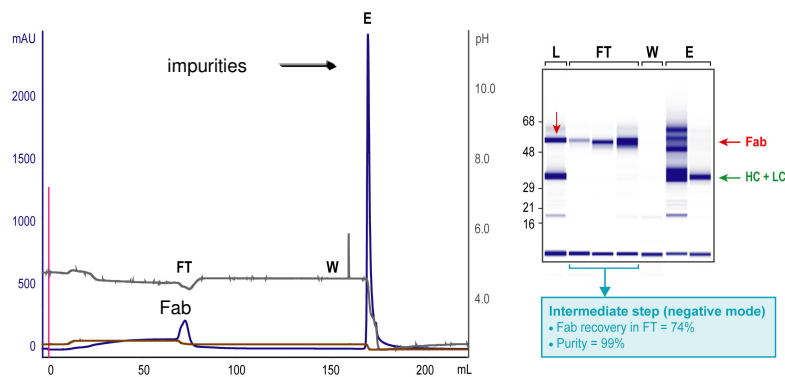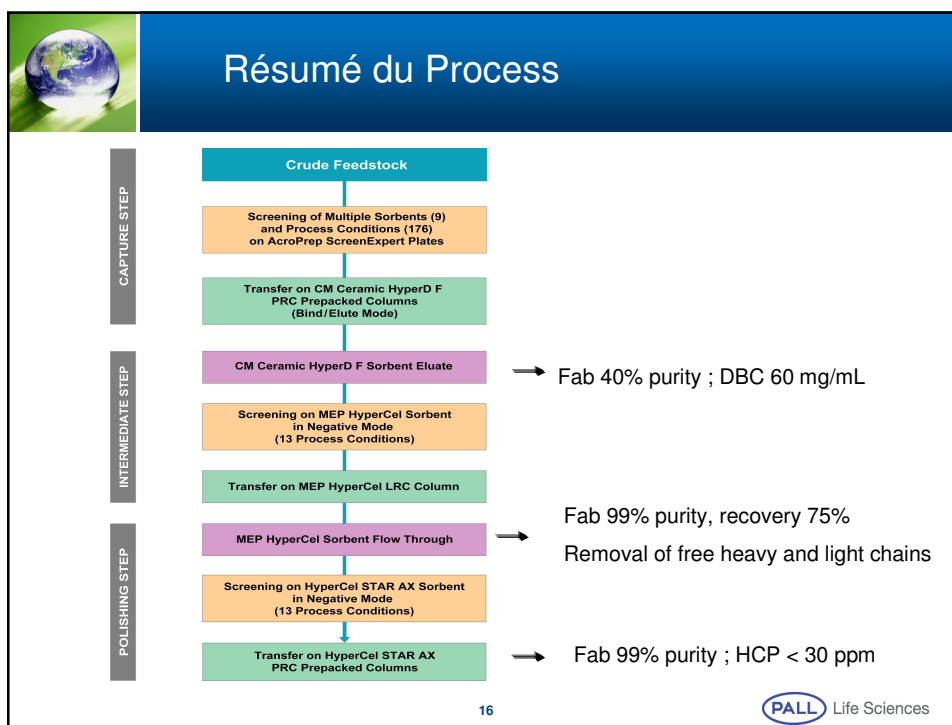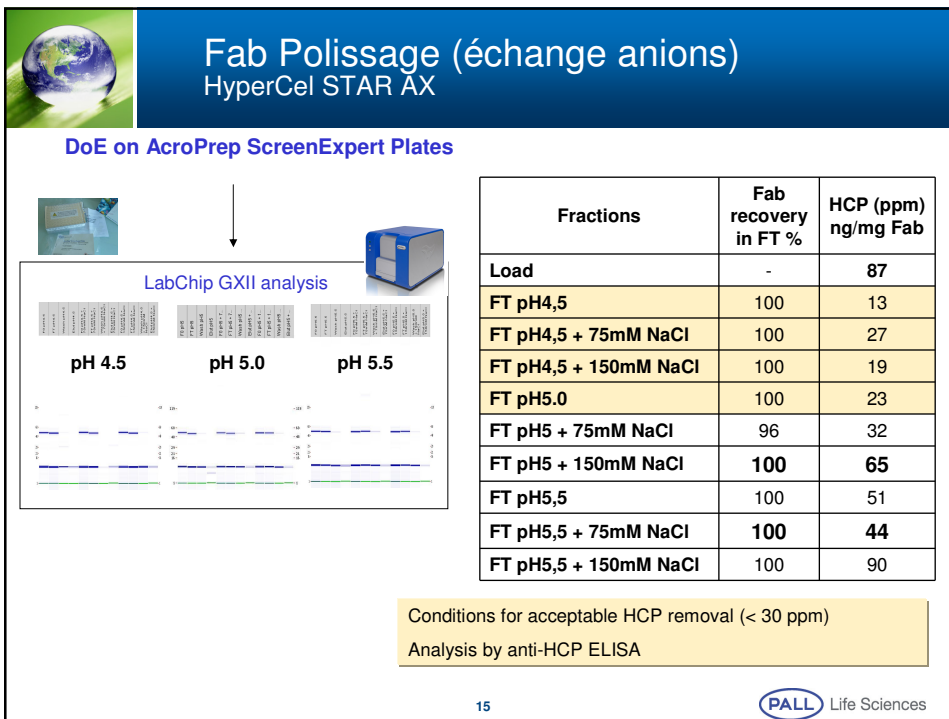Equilibration: 50 mM Na acetate pH 5.2
Loading: 100 mL feedstock 1/10 at pH 5.2
Washing: 50 mM Na acetate pH 5.2
Elution: 50 mM Na acetate pH 3.0
Flow rate: 150 cm/h – Residence Time = 5 min

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Publications

- [1] High throughput screening of mixed-mode sorbents and optimization using pre-packed lab-scale columns for the purification of the recombinant allergen rBet v 1a. V Brenac Brochier *et al.*, **J. Chromatogr. B** (2009).
- [2] Rapid screening of purification strategies for the capture of a human recombinant F(ab')₂ expressed in baculovirus-infected cells using a micro-plate approach and SELDI-MS. J Pezzini *et al.*, **J. Chromatogr. B** (2009).
- [3] Expression optimization and purification process development of an engineered soluble recombinant mouse linker of activation of T cells using SELDI-MS. V Brenac *et al.*, **Prot. Express. and Purif.** (2006).

**ENABLING A
GREENER
FUTURESM**

www.pall.com/green

This presentation is the copyright work product of Pall Corporation and no portion of this presentation may be copied, published, performed, or redistributed without the express written authority of a Pall corporate officer.

© 2013 Pall Corporation